

APEX MICROSCOPES – 120 (AM-120)

60-120x LED Lighted Pocket Microscope

User and Safety Manual

Congratulations on selecting your new LED Lighted Pocket Microscope!
In order to achieve optimum performance, please follow instructions for proper use and care.

PRODUCT SPECIFICATION


Magnification	60x – 120x
Optical Zoom	2x
Exit Pupil Diameter	10 mm – 14 mm
Illumination	LED
Battery	1 × AA
Weight	0.13 lbs
Dimensions	2.3" × 1.2" × 3.7"

WARNING

- To avoid permanent eye damage:**
 - Do not look at the sun with the microscope
 - Do not place the illuminated microscope directly on the human eye
- Remove the battery before storing the microscope for extended periods
- The device may become warm after extended use. This is normal and should not be considered a defect.

INSTALLING THE BATTERY

- Slide battery cover to open.
- Insert 1 AA alkaline battery and slide back battery cover.

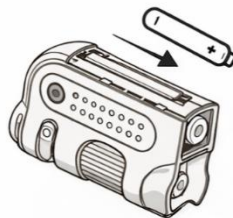

⚠ WARNING

- Use only alkaline battery.
- Keep away from fire or excessive heat.
- Inspect battery for corrosion or leakage.
- Insert in direction indicated.
- Rechargeable battery not recommended.

USING THE MICROSCOPE

- Place viewing object directly below Objective Lens.
- Press LED switch to turn light On and Off.
- Set Zooming Dial to lowest magnification and rotate Focusing Ring slowly until image is clear. Adjust zoom and refocus as needed.



Note: To work properly, microscope must be placed flush to viewing object.

CUSTOMER SUPPORT

For technical assistance or product support, please contact our team. We are committed to providing prompt and reliable service.

✉ info@eu.apexinternational.com

🌐 www.apexinternational.com